



DATA SHEET

No Restriction

Title:	Afinion™ Data Connectivity Converter Interface HL7		
Version no.:	3.0	Pages:	5
Main project:	Afinion™ Data Connectivity Converter		
Sub project:	Data Sheets	Doc. ID:	TP09/40/602/609/ 682-4486
Prepared by:	Vegar Kjekshus	Date:	30.06.2010
Approved by:	Marianne Solberg (SOLB) 02.07.2010 06:36:06		

1. INTRODUCTION

This document describes the Data Connectivity Protocol (patient and control records) for LIMS connectivity based on the HL7 (health level 7) protocol in the ADCC (Afinion™ Data connectivity converter) for the Afinion™ AS100 Analyzer. It describes the protocol and the format of the records returned from the Analyzer. It also gives examples and highlights issues to be especially addressed; all needed by the programmer that shall interface to this protocol on the LIMS side.

1.1 Revision

Revision	Date	Author	Comment
1	28.01.2010	DZe/ TUh/ ESc	First version
2	10.06.2010	VEKJ	Ch. 1.1 Deleted unused cells. Ch. 4.2, 4.5, 4.6,4.7,4.8, 5.1 Corrected and changed examples Ch. 5.1 Added example.
3	30.06.2010	VEKJ	Ch.4.8, 5.1.3 Updated with precaution and example for ACR out of range results

2 PHYSICAL TRANSMISSION OF MESSAGES

2.1 HL7 Socket transfer

Each HL7 message will be transmitted as follows:

Sender (PCC)	Direction	Receiver (LIS)
<VT>HL7-ORU-O01-message<FS><CR>	-->	
	<--	<VT>HL7-ACK-message<FS><CR>

wherein:

<VT> ... ASCII 0x0B
 <FS> ... ASCII 0x1C
 <CR> ... ASCII 0x0D

3 MESSAGE STRUCTURE

3.1 HL7 message structure

The HL7 high level message structure is based on Version 2.4 of the Health Level Seven (HL7) Standard for electronic data exchange in all healthcare environments.

3.2 Patient measurement results to LIS

The Afinion will transmit the following events for patient measurement results:

Message type	Event code	Description
ORU	R01	Unsolicited transfer of patient results

Message structure:

HL7 segment	Description
MSH	Message Header
PID	Patient Identification
PV1	Patient Visit
OBR	Observation Request
OBX	Observation Result

4 SEGMENTS

4.1 Legend

- Field name: according to Health Level Seven, version 2.4
- Interpretation: additional description
- Req.: F ... fix value, C ... configured value via web interface, A ... data comes from analyzer, X ... calculated values (e.g. date/time), R ... required from LIS, O ... optional from LIS
- HL7 Field: Nr of HL7 field
- Source of data: Field of the Afinion data record, where the data comes from.

4.2 MSH (H)

Field name	Interpretation	Req.	HL7 Field	contents
Field separator	Field separator byte	F	1	
Encoding characters	Other field separator characters	F	2	
Sending application	Model name (always "Afinion AS100")	F	3	
Receiving application	Name of the receiving application / dept. (configurable)	C	5	Configured value
Receiving facility	Receiving process / institution within the dept. (configurable)	C	6	Configured value
Date / time of message	date and time of message creation	X	7	current time stamp
Message type	always "ORU"	F	9.1	
Event type	always "R01"	F	9.2	
Message Control ID	Consecutive number starting with 1000	X	10	1000
Processing ID	P ... patient measurement results Q ... quality control results	A	11	P for record,patient@ Q for record,control@
Version ID	HL7-version used	F	12	HL7: "2.4"
Accept Acknowledgement Type	always "AL"	F	15	AL
Application Acknowledgement Type	always "NE"	F	16	NE
Character Set	always "8859/1"	F	18	8859/1

HL7-Example: MSH|^~\&|Afinion AS100||EPR|KH-1|20100610131643||ORU^R01|1048|P|2.4|||AL|NE||8859/1

4.3 EVN

EVN is not supported for patient measurement result export to LIS.

4.4 MSA

Field name	Interpretation	Req.	HL7 Field	Contents
Acknowledgement code	AA, CA will be accepted as acknowledgement from LIS	R	1	AA
Message Control ID	Verification of message control ID will not be performed.	O	2	
Text Message	Error text will not be analyzed.	O	3	
ErrorCondition	Error code, if Acknowledgement code is neither AA nor CA, Error condition will be saved into the log memory.	O	6	

HL7-Example: MSA|AA|117715205|||F|

4.5 PID (P)

Field name	Interpretation	Req.	HL7 Field	contents
Set ID - Patient ID	PID segment number	F	1	
Patient Identifier List	(local) patient ID	A	3	P- ID of header (configurable, to use PID or visitnumber)

HL7-Example: PID|1||43|

5 EXAMPLES

5.1 ORU-Message (HL7)

5.1.1 Example 1

```
MSH|^~\&|Afinion AS100||EPR|KH-1|20100610131643||ORU^R01|1048|P|2.4|||AL|NE||8859/1
PID|1|||
PV1|1|||43|
OBR|1||3|CRP|||N|||
ORH|||^^10124809|F|
OBX|1|NM|CRP||16|mg/L|||F|||AS0007962|20100608142352|
```

5.1.2 Example 2

```
MSH|^~\&|Afinion AS100||EPR||20100608185608||ORU^R01|1006|P|2.4|||AL|NE||8859/1
PID|1||55|
PV1|1|||
OBR|1||1|ACR|||N|||ORH|||^^10142193|F|
OBX|1|NM|ACR||0.5|mg/g|||F|||AS0007962|20100608140517|
OBX|2|NM|Alb||8.0|mg/L|||F|||AS0007962|20100608140517|
OBX|3|NM|Creat||17.4|mg/dL|||F|||AS0007962|20100608140517|
```

5.1.3 Example 3

This example shows an ACR measurement where the ACR result must be interpreted as **not valid** even if abnormal flags are missing since the albumin (Alb) result is below measurement range.

```
MSH|^~\&|Afinion AS100||EPR||20100616123242||ORU^R01|1008|P|2.4|||AL|NE||8859/1
PID|1||2|
PV1|1|||
OBR|1||2|ACR|||N|||ORH|||^^10142193|F|
OBX|1|NM|ACR||5.6|mg/g|||F|||AS0007962|20100608140536|
OBX|2|NM|Alb||4.1|mg/L||<||F|||AS0007962|20100608140536|
OBX|3|NM|Creat||33.0|mg/dL|||F|||AS0007962|20100608140536|
```